High sensitivity type (integrated with back-thinned type CCD image sensor)

TG series mini-spectrometers are polychromators integrated with optical elements, an image sensor and a driver circuit. Light to be measured is guided into the entrance port of TM series through an optical fiber and the spectrum measured with the built-in image sensor is output from the USB port to a PC for data acquisition. C9404CA, C9404CAH and C9405CA are high sensitivity mini-spectrometers employing a back-thinned type CCD image sensor. Their sensitivity is about two orders of magnitude higher than CMOS types making them even more ideal for low-light-level measurement. C9404CAH is high resolution type (resolution: 1 nm Typ.). Their products come supplied with free sample software that allows setting measurement conditions, acquiring and saving data, and displaying graphs. Driver software and DLL are also supplied as accessory items to allow the users to configure their own measurement software.

#### **Features**

- Integrated with back-thinned type CCD image sensor: Sensitivity is about two orders of magnitude higher than CMOS types.
- High resolution 1 nm (C9404CAH)
- High throughput due to transmission grating made of quartz
- Highly accurate optical characteristics
- Easy to install into equipment
- Wavelength conversion factor is recorded in internal memory \*1

#### Optical characteristics

Optical characteristics					
Parameter	TG-UV-CCD		TG-SWNIR- CCD	Unit	
	C9404CA	C9404CAH	C9405CA		
Spectral response range	200 to 400		500 to 1100	n m	
Spectral resolution Max.	3	1 *3	5 *4	n m	
(Spectral response half width) *2	3		3	11111	
Wavelength reproducibility *5	±0.1		±0.2	n m	
Wavelength temperature	0.02			nm/°C	
dependence	0.02			IIIII/ C	
Spectral stray light *2, *6	-35			dB	

#### ■ Electrical characteristics

Parameter	Specification	Unit
A/D conversion	16	bit
Integration time	10 to 10000	m s
Interface	USB1.1	-
USB bus power current consumption	100	m A
External power supply	5	V

#### ■ General ratings / Absolute maximum ratings

Specification			Unit
125.7 (W) × 115.7 (D) × 75 (H)			m m
Back-thinned type CCD image sensor (S10420-1006)			-
1024			pixels
140 × 500	10 × 1000	70 × 800	μm
0.11 0.22		0.22	-
SMA905D			-
+5 to +40			°C
-20 to +70			°C
	125.7 (W Back-thingsens 140 × 500	125.7 (W) × 115.7 (D)  Back-thinned type CC sensor (S10420-1 1024 140 × 500   10 × 1000 0.11 SMA905D +5 to +40	125.7 (W) × 115.7 (D) × 75 (H)  Back-thinned type CCD image sensor (S10420-1006)  1024  140 × 500

- \*1: A conversion factor for converting the image sensor pixel number into a wavelength is recorded in the module. A calculation factor for converting the A/D converted count into the input light intensity is not provided.
- \*2: Depends on the slit opening. Values were measured with the slit listed in the table "


  General ratings /Absolute maximum ratings".
- \*3: Typical
- \*4: λ=550 to 900 nm
- \*5: Measured under constant light input conditions
- \*6: When monochromatic light of the following wavelengths is input, spectral stray light is defined as the ratio of the count measured at the input wavelength, to the count measured in a region of the input wavelength ±20 nm (C9404CA, C9404CAH) or ±40 nm (C9405CA).
  - C9404CA, C9404CAH: 300 nm, C9405CA: 800 nm
- \*7: Entrance slit aperture size
- \*8: No condensation

#### PRELIMINARY DATA Jan. 2007

# **Applications**

- Low-light-level measurement such as fluorescence measurement
- Detection of saccharic acids in foods
- Taste analyzers
- Evaluation of light source characteristics such as UV light source


### Comparison of CCD type and CMOS type

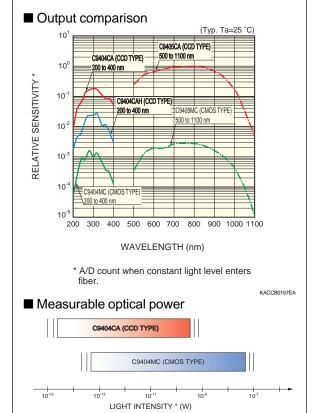

\* Intensity of light incident on mini-spectrometer through slit is constant. (λ=300 nm, Integration time 10 to 10000 ms)




# ■ Spectral resolution vs. wavelength

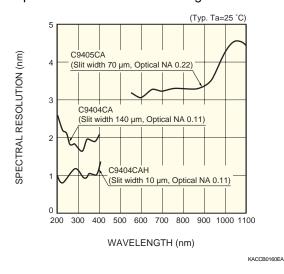


# ■ Optical component layout

TM series mini-spectrometers use a transmission holographic grating made of quartz and precision optical components arranged on a rugged optical base, making it possible to deliver high throughput and highly accurate optical characteristics.

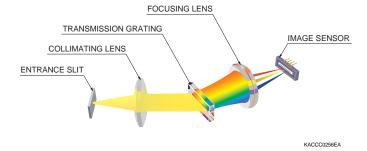


■ Connection example (transmission light measurement)
Light to be measured is guided into the entrance port of TG
series through an optical fiber and the spectrum measured
with the built-in image sensor is output through the USB port
to a PC for data acquisition. There are no moving parts inside
the unit so stable measurements are obtained at all times. An
optical fiber that guides light input from external sources allows a flexible measurement setup.

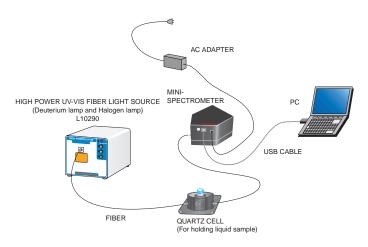

KACCC0355EA

## ■ Measurement example

Line spectra from low-pressure mercuy lamp were measured with C9404CA (TG-UV-CCD), C9404CAH (TG-UV-CCD).

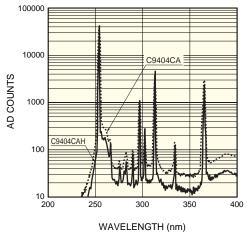

KACCB0166EA

## ■ Dedicated software package (supplied with unit)

Installing the dedicated software package (containing sample software, device driver, DLL)\*9 into your PC allows running the following basic tasks:

- · Measurement data acquisition and save
- · Measurement condition setup
- Module information acquisition (wavelength conversion factor, polychromator type, etc.)
- · Graphic display
- · Arithmetic operation

Pixel number to wavelength conversion

Dark subtraction

Comparison calculation with reference data


(transmittance, reflectance)

Gaussian approximation

(peak position and count, FWHM)

Note: Two or more mini-spectrometers can be connected and used with one PC simultaneously.

\*9: Compatible OS: Microsoft Windows Professional Edition 2000 (SP3 or later) and XP (SP1a or later)

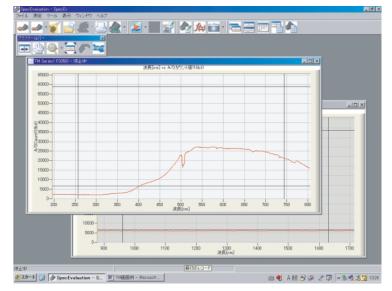


Device driver and DLL for controlling hardware are also provided.

You can develop your own measurement programs by using a software development environment that includes Microsoft Visual C++ and Visual Basic.\*<sup>10</sup> The DLL provides functions such as USB port open/close, measurement condition setup, measurement data and module information acquisition.

\*10: Operation of the device driver and DLL has been verified only with Microsoft Visual C++® and Visual Basic®.

Microsoft Visual C++ and Microsoft Visual Basic are either registerd trademarks or trademarks of Microsoft Corporation in the United States and other countries.

#### ■ Dimensional outline (unit: mm)

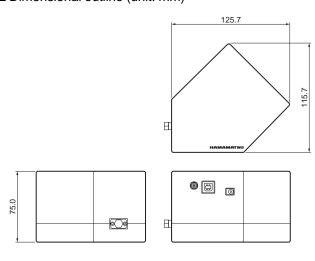

KACCA0202EA


#### Accessories

- · USB cable
- · Dedicated software (sample software, device driver, DLL)
- · AC adapter (for power supply)

## ■ Options (sold separately)

Optical fibers for light input

Type. No	Product name	Applicable mini-spectrometer	Core diameter (µm)	Specification
A9762-01	Fiber for UV/visible range (resistance to UV)	C9404CA (TG-UV-CCD) C9404CAH (TG-UV-CCD)	600	N.A.=0.22, length 1.5 m, connectorized SMA905D at both ends
A9763-01	Fiber for visible/near infrared range	C9405CA (TG-SWNIR-CCD)	600	

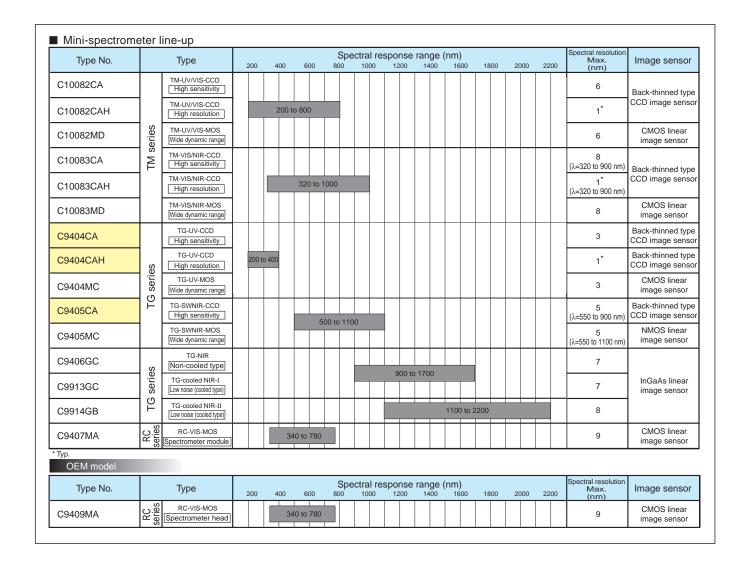

HAMAMATSU Information Specificat

Information furnished by HAMAMATSU is believed to be reliable. However, no responsibility is assumed for possible inaccuracies or omissions.

Specifications are subject to change without notice. No patent rights are granted to any of the circuits described herein. ©2007 Hamamatsu Photonics K.K.

HAMAMATSU PHOTONICS K.K., Solid State Division

1126-1 Ichino-cho, Higashi-ku, Hamamatsu City, 435-8558 Japan, Telephone: (81) 53-434-3311, Fax: (81) 53-434-5184, www.hamamatsu.com U.S.A.: Hamamatsu Corporation: 360 Foothill Road, P.O.Box 6910, Bridgewater, N.J. 08807-0910, U.S.A., Telephone: (1) 908-231-0960, Fax: (1) 908-231-1218 Germany: Hamamatsu Photonics Deutschland GmbH: Arzbergerstr. 10, D-82211 Herrsching am Ammersee, Germany, Telephone: (49) 08152-3750, Fax: (49) 08152-2658

France: Hamamatsu Photonics France S.A.R.L.: 19, Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: 33-(1) 69 53 71 00, Fax: 33-(1) 69 53 71 10
United Kingdom: Hamamatsu Photonics UK Limited: 2 Howard Court, 10 Tewin Road, Welvyn Garden City, Hertfordshire AL7 18W, United Kingdom, Telephone: (44) 1707-294888, Fax: (44) 1707-325777
North Europe, Hamamatsu Photonics Wick Sprides/Sept. 12 SE-1374 (1 Solns, Sweden, Telephone; (46) 8-500-031-01

North Europe: Harnamatsu Photonics Norden AB: Smidesvägen 12, SE-171 41 Solna, Sweden, Telephone: (46) 8-509-031-00, Fax: (46) 8-509-031-01 Italy: Harnamatsu Photonics Italia S.R.L.: Strada della Moia, 1/E, 20020 Arese, (Milano), Italy, Telephone: (39) 02-935-81-733, Fax: (39) 02-935-81-741